

Washing Machine for Bottles WT-2000




WT turn table type cartridge bathing device

Highly reliable cleansing, cartridge supply by turn table, various size of cartridge and wide selection of production range

WT device is mainly used for cleaning that requires no ultrasonic cleaning. Container is supplied through revolving table with the size of 1,000mm maximum diameter. The container is arranged in a row and transported to moving gripper through Inlet star wheel. Then, moves to the next process as it is outputted through outlet star - wheel.

Processing sequences:

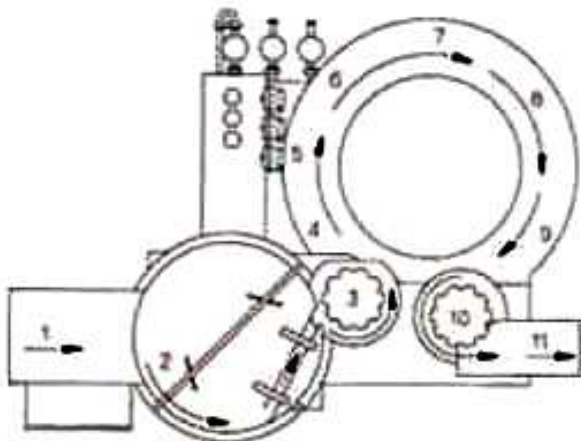

1. Supply table
2. Revolving table
3. Inlet Star Wheel
4. Catch the container then turn 180
5. Clean the internal/external parts of the container with distilled water
6. Clean the inside of the container with compressed air
7. Clean the inside of the container with distilled water
8. Clean the internal/external parts of the container with compressed water
9. Silicon treatment to the container (option)
10. Outlet star Wheel
11. Output to next processing stage



Container supply



Send container and cleaning



Container output



Air & cleaning water filter



Input gauge & control valve



Distilled water supply tank and pump

Technical Data:

Description		MODEL WT
		2000
Number of Gripper		40
Output (Vials/min : 10mℓ)		200
Electricity	Installed Power	220/380V, 50/60Hz. 16KW
Consumption	Main Drive	0.75KW
	Recycling Pump	2.2 KW
	Boost Pump	1.5 KW
	Heater (option)	9 KW
	Ultrasonic Generator	None
Dimension	Length	2700 mm
	Width	1980 mm
	Hight	1380 mm